



Mechanical Pipettes

MicroPette plus

121°C
FULLY
Autoclave

Features

- Fully autoclavable
- Ergonomic design provides excellent operating experience
- Easy-to-read volume display
- The pipettes cover the volume range from 0.1µL to 10mL
- Easy calibration and maintenance
- Manufactured from innovative materials
- Each MicroPette Plus supplied with an individual calibration certificate according to ISO8655



MicroPette plus multi-channel

121°C
FULLY
Autoclave

Features

- 8 and 12 channel pipettes are appropriate for 96-well plates
- Dispensing head rotates for effortless pipetting convenience
- Individual piston and tip cone assemblies allows for easy repair and maintenance
- Compound material-made tip cone secures high sealing performance
- Compatible with most universal tip brands

Push button for volume adjustment

Comfortable Finger support

Low force tip ejector for easy operation

Spring loaded tip cones for easy cleaning and maintenance

High chemical-resistance materials to secure long lifetime

8 channels

Fully autoclavable

The MicroPette Plus pipettes are fully Autoclavable enable easy cleaning to reduce the risk of contamination. Steam autoclaving can be performed at 121 °C, 1 bar for 20 minutes. After autoclaving the pipette must be cooled down and left to dry for 12 hours before use.

It is recommended to check the performance of the pipette after each autoclaving. Greasing and sealing of the pipette piston after every 10th autoclaving will secure an enhanced function.

