
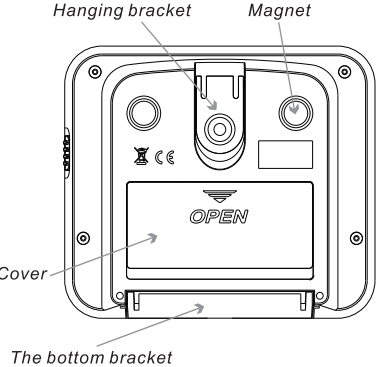


User's Guide

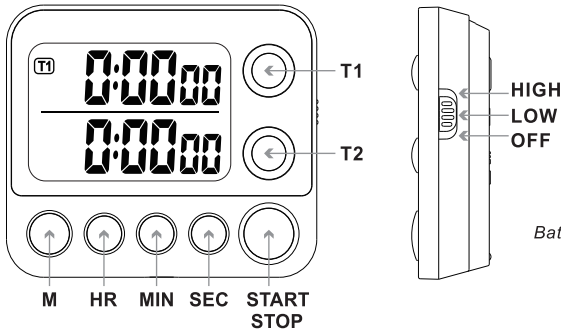

- T1** button ----- To switch to control the first set timer T1.
- T2** button ---- To switch to the control of the second set timer T2.
- M** button ----- To set the memory time, return to the memory time.
- HR** button ----- To adjust "hour" value.
- MIN** button ----- To adjust "minute" value.
- SEC** button ---- To adjust "second" value.
- START / STOP** button ---- To start or pause the timer.

Set memory time:

You can quickly re-use a time setting as memory time.

Set as follows:

- 1 When the timer is paused, press and hold the **M** button until the value flashes.
 - 2 Press the **HR** button to increase "hour" value, press and hold to increase quickly.
 - 3 Press the **MIN** button to increase "minute" value, press and hold to increase quickly.
 - 4 Press the **SEC** button to increase "second" value, press and hold to increase quickly.
- * Press the **HR** and **MIN** buttons to clear the timer.
 - 5 Press the **M** button to finish the default time setting.
 - * Whenever you want to use the memory time, press the **M** button to bring up the memory time you set, when the timer is paused.

Use the stopwatch - >>

- * Press the **T1** or **T2** button to switch to the **T1** or **T2** timer you want to use.
- 1 Press the **HR** and **MIN** buttons to clear the timer when the timer is paused.
 - 2 Press the **START / STOP** button to start the stopwatch immediately.
- * The timer is accurate to centisecond within one hour, and accurate to seconds after one hour. The timer counts up to 99 hours 59 minutes 59 seconds.

Control alarm - >>

The fluctuation switch can be switched to corresponding position to adjust the volume level of the alarm.

HIGH: High volume **LOW:** low volume **OFF:** Alarm off


How to place the product - >>

- * You can use the hanging bracket to hang it on the wall or other places.
- * The product can stand upright on the table with the bottom bracket unfolded.
- * This product comes with a magnet, it can be attached onto the metal surface of the iron box or refrigerator.

How to replace the battery - >>

When the power is low, the display becomes dim or unclear. Replace the battery when the battery power is low.

- 1 Hold and push the battery cover towards the direction indicated by "▲" to open the battery cover.
- 2 Remove the old battery, and put on a new battery according to the correct polarity direction.
- 3 Replace the battery cover.

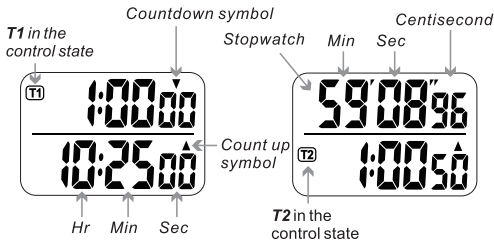

Timer Instructions - >>

* Press the **T1** or **T2** button to switch to **T1** or **T2** timer you want to use.

Set the countdown time:

- 1 Press the **HR** button to increase "hour" value, press and hold to increase quickly.
 - 2 Press the **MIN** button to increase "minute" value, press and hold to increase quickly.
 - 3 Press the **SEC** button to increase "second" value, press and hold to increase quickly.
 - 4 Press the **HR** and **MIN** buttons to clear the timer.
- * You can set up to 99 hours 59 minutes 59 seconds of the countdown time.

Start the countdown

* After setting the countdown time, press the **START / STOP** button to start the countdown, with the "▼" flashing; Press the **START / STOP** button again to pause countdown, with the "▼" stop flashing; Press the **START / STOP** button again to continue the countdown.

* When the countdown is in progress, the corresponding button lamp (green) of the timer you are using blinks.

After the end of the countdown

- * When the timer counts down to zero, the count-up symbol "▲" flashes, the timer will count up from zero, and the alarm will sound for a minute (if the alarm is turned on using fluctuation switch), and the corresponding indicating lamp (red) will flash for a minute.
- * Press the **START / STOP** button to return immediately to the set countdown time.